



INSTALLATION, OPERATION & MAINTENANCE INSTRUCTIONS

1455 Kleppe Lane ♦ Sparks, NV 89431-6467 ♦ (775) 359-4712 ♦ Fax (775) 359-7424
E-mail: haws@hawsco.com ♦ website: www.hawsco.com

No. 0510002288 (2)

Model 9006 PORTABLE EYEWASH STAND

NOTE TO INSTALLER: Please leave this information with the Maintenance Department.

LIMITED WARRANTY

HAWS warrants that all of its products are guaranteed against defective material or poor workmanship for a period of **one year from the date of shipment**. The foregoing notwithstanding, *HAWS* warrants certain specific products or components thereof for an adjusted period. A list of these excepted products and components and a description of their respective warranty terms may be found here: www.hawsco.com/warranty. *HAWS'* liability under this warranty shall be discharged by furnishing without charge F.O.B. *HAWS* factory any goods, or part thereof, which shall appear to the Company upon inspection to be of defective material or not of first-class workmanship, provided that a claim is made in writing to Company within a reasonable period after receipt of the product. Where claims for defects are made, the defective part or parts shall be delivered to the Company, prepaid, for inspection. *HAWS* shall not be liable for the cost of repairs, alterations or replacements, or the labor required to implement them, or for any expense connected therewith made by the owner or owner's agents, except upon written authority from *HAWS*. *HAWS* shall not be liable for any damages caused by defective materials or poor workmanship, except for replacements, as provided above. Buyer agrees that *HAWS* has made no other warranties either express or implied in addition to those above stated, except that of title with respect to any of the products or equipment sold hereunder and that *HAWS* shall not be liable for general, special, incidental, or consequential damages claimed to arise under the contract of sale.

Equipment manufactured by *HAWS* is warranted to function if installation and maintenance instructions provided are adhered to. The units also must be used for the purpose for which they were intended. Any *HAWS* emergency equipment is intended to supplement first-aid treatment. Due to widely varying conditions, *HAWS* cannot guarantee that the use of this emergency equipment will prevent serious injury or the aggravation of existing or prior injuries.

EXCEPT AS EXPRESSLY STATED HEREIN, HAWS HEREBY DISCLAIMS ALL WARRANTIES, WHETHER EXPRESS OR IMPLIED, ARISING BY LAW OR OTHERWISE, INCLUDING WITHOUT LIMITATION, ANY IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. THIS WARRANTY MAY NOT BE MODIFIED OR EXTENDED WITHOUT THE WRITTEN CONSENT OF HAWS. REMEDIES AND REPLACEMENTS STATED HEREIN ARE EXCLUSIVE. IN NO EVENT SHALL HAWS BE LIABLE FOR ANY SPECIAL, PUNITIVE, CONSEQUENTIAL OR INCIDENTAL DAMAGES TO ANY PERSON INCLUDING BUT NOT LIMITED TO DAMAGES FOR LOSS OF USE OR PROFITS, SUBSTITUTE PRODUCTS OR COSTS, PROPERTY DAMAGE, OR OTHER MONETARY LOSS.

SHOULD YOU EXPERIENCE DIFFICULTY WITH THE INSTALLATION OF THIS MODEL PLEASE CALL:

TECHNICAL SUPPORT: 1-800-766-5612
FOR CUSTOMER SERVICE: 1-888-640-4297

LOCATION OF UNIT: The Model 9006 is intended for use indoors on smooth, level ground. Refer to your eyewash installation instructions for additional considerations when locating the unit.

COMPATIBLE EYEWASH PRODUCTS: The Model 9006 is designed to support an eyewash with a maximum weight of 175lb and is compatible with a wide range of portable eyewash models. Verify your eyewash mounting requirements and refer to the dimensions provided in the Model 9006 installation instructions to determine compatibility.

INSTALLATION PROCEDURE

GENERAL NOTES:




- 1. REFER TO INSTALLATION INSTRUCTIONS ON THE FOLLOWING PAGES.**
- 2. AFTER INSTALLATION, CHECK EYEWASH FUNCTION TO ENSURE THAT THE UNIT IS STILL IN COMPLIANCE WITH THE REQUIREMENTS OF PERTINENT CODES AND STANDARDS (E.G., DIMENSIONS, ACTIVATION, FLOW CHARACTERISTICS, ETC.).**
- 3. ALWAYS EMPTY AND/OR REMOVE AN INSTALLED EYEWASH PRIOR TO RELOCATING THE MODEL 9006 TO AVOID DAMAGE OR INJURY DUE TO TIPPING.**


*For more information about Haws products, see our website: www.hawsco.com



NOTES:

1. HOLD ROUGH-IN DIMENSIONS $\pm 1/4"$ (6.4mm).
2. WHEN INSTALLING THIS UNIT, LOCAL, STATE, OR FEDERAL CODES SHOULD BE ADHERED TO. FOR INSTALLATION HEIGHTS OTHER THAN SHOWN, DIMENSIONS MARKED (*) MUST BE ADJUSTED ACCORDINGLY. INSTALL A TRAP IN THE WASTE LINE AND A SCREWDRIVER STOP IN THE SUPPLY LINE IF REQUIRED.
3. TO COMPLY WITH ANSI Z358.1-2014 FOR EMERGENCY EYEWASH AND SHOWER EQUIPMENT:
 - UNIT SHALL BE INSTALLED IN ACCORDANCE WITH MANUFACTURER'S INSTRUCTIONS AND ACCEPTABLE PLUMBING PRACTICES.
 - EYEWASH STREAMS SHALL BE POSITIONED SUCH THAT STREAMS ARE 15.3 CM (6 IN.) MINIMUM FROM THE WALL OR NEAREST OBSTRUCTION AND MEET ANSI GAUGE REQUIREMENTS BETWEEN 83.8 & 134.6 CM (33-53 IN) ABOVE FLOOR.

(NOTES CONTINUED ON NEXT SHEET)

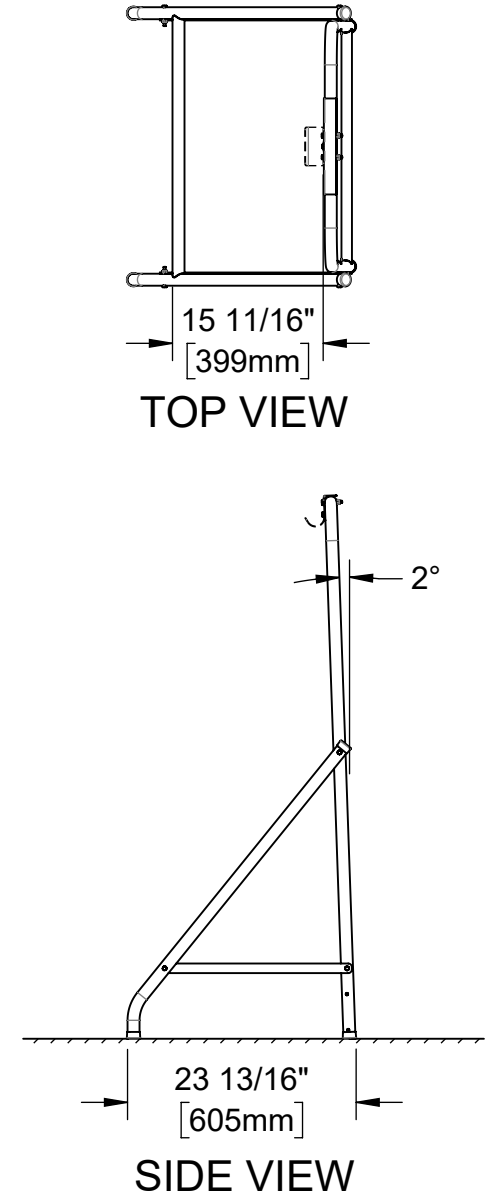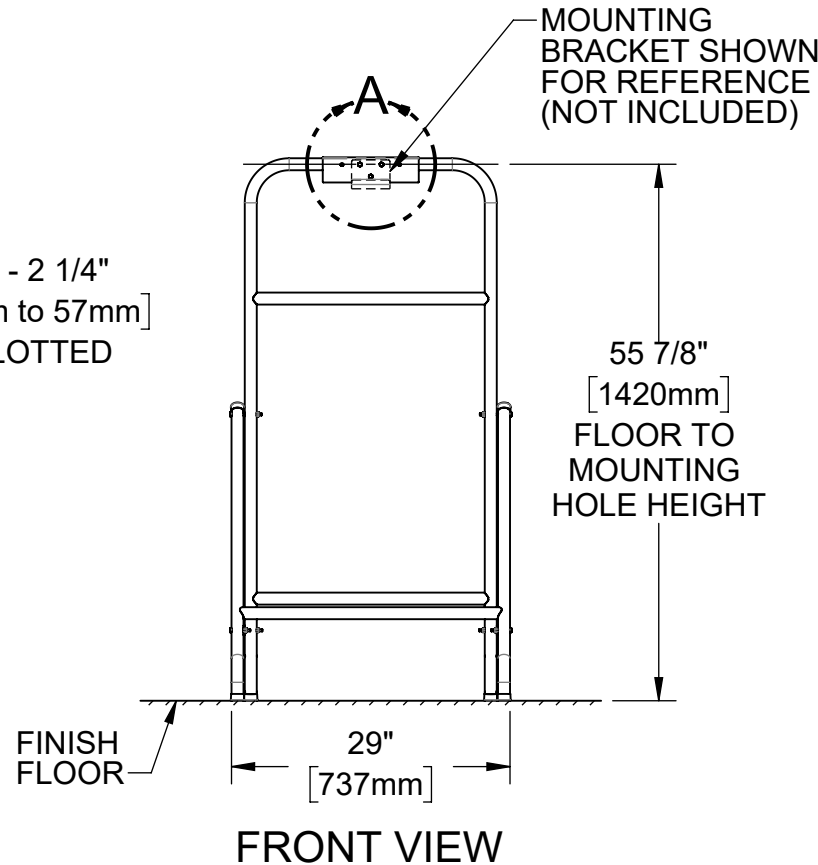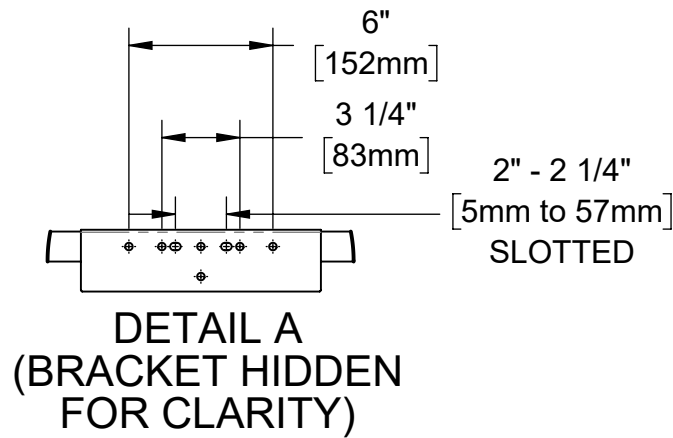

				1455 KLEPPE LANE SPARKS, NEVADA 89431 (775) 359-4712 FAX (775) 359-7424 E-MAIL: HAWS@HAWSO.COM WEBSITE: WWW.HAWSO.COM			
ECN: 6210	REV. ECN: 6305	BY: JU	MODEL(S)	9006, PORTABLE EYEWASH STAND			PART NUMBER
DRAWN: JU	DATE: 01/13/25	CHKD: JL					0510002288.D
APPROVED: JU	DATE: 07/29/25						REVISION
							2
SCALE: 1:20				DRAWING TYPE: INSTALLATION		SIZE: A	SHEET 1 OF 3

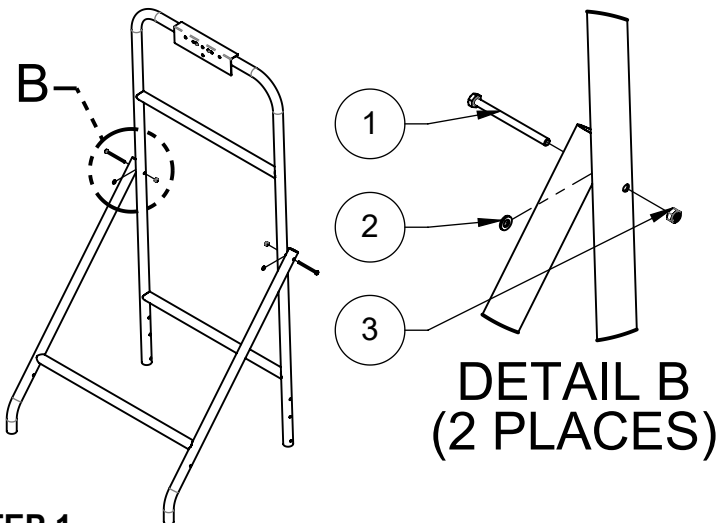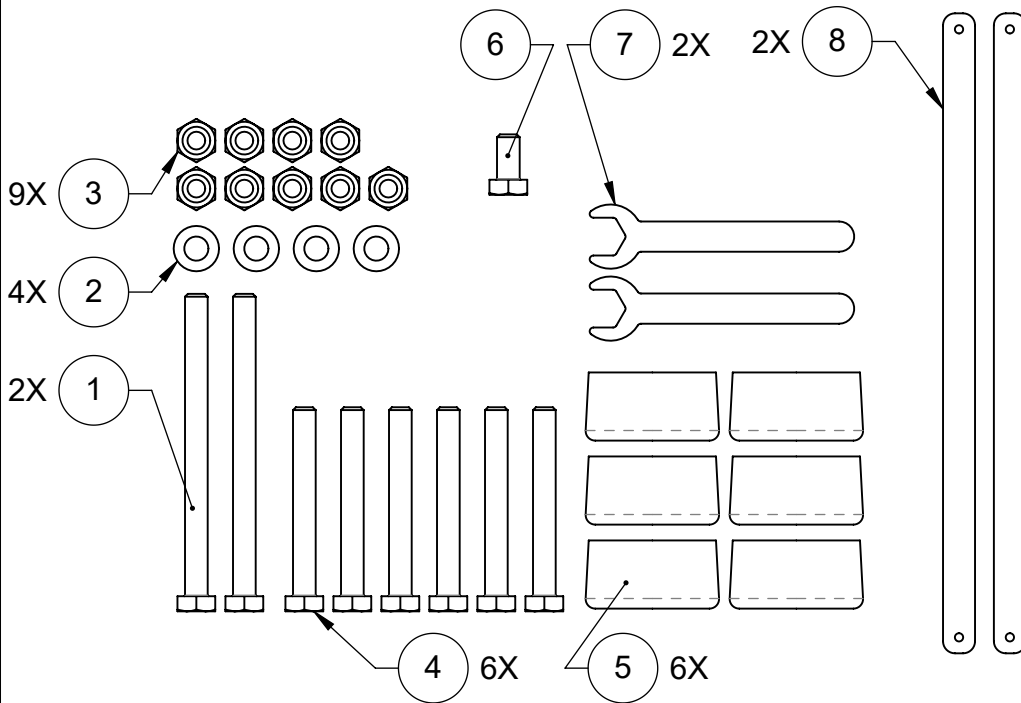


**DETAIL B
(2 PLACES)**

STEP 1
CONNECT THE TUBULAR FRAMES USING LONG SCREWS (ITEM 1), WASHERS (ITEM 2), AND LOCK NUTS (ITEM 3). WASHERS (ITEM 2) ARE INSTALLED BETWEEN THE FRAME PIECES.

8	BRACKET, LOWER BRACE	2
7	WRENCH, OPEN END, 10MM	2
6	SCREW, M6 X 12MM, HEX	1
5	CAP, SLIP ON, PLASTIC	6
4	SCREW, M6 X 50MM, HEX	6
3	NUT, M6, W/ NYLON INSERT	9
2	WASHER, M6	4
1	SCREW, M6 X 80MM, HEX	2
ITEM	DESCRIPTION	QTY
INCLUDED HARDWARE		

NOTES CONTINUED:

- TORQUE ALL NUTS AND BOLTS TO 26.5 LB-IN [3 Nm]. OVER-TORQUEING HARDWARE MAY DAMAGE STEEL TUBING.
- RECOMMENDED TOOLS:
 - 10MM SOCKET AND SOCKET DRIVER.

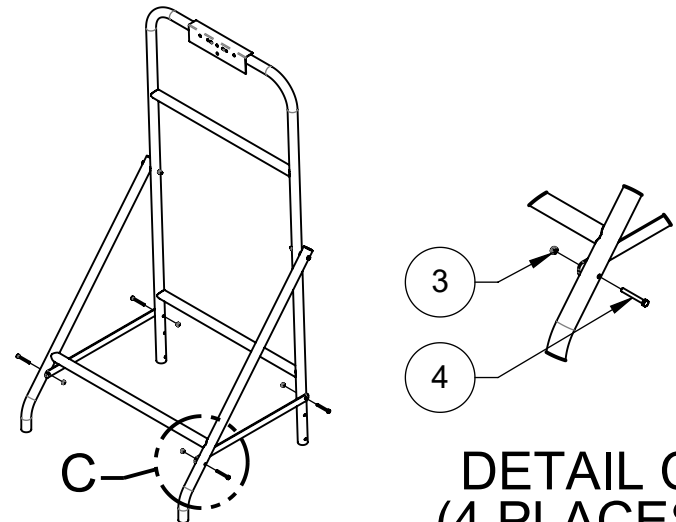


**DETAIL C
(4 PLACES)**

STEP 2
INSTALL THE TWO FLAT BRACKETS (ITEM 8) AS SHOWN, USING MEDIUM SCREWS (ITEM 4) AND LOCK NUTS (ITEM 3). NOTE THESE BRACKETS ATTACH ON THE INSIDE SURFACE OF THE FRONT FRAME AND OUTER SURFACE OF THE VERTICAL FRAME.




1455 KLEPPE LANE
SPARKS, NEVADA 89431
(775) 359-4712 FAX (775) 359-7424
E-MAIL: HAWS@HAWSO.COM
WEBSITE: WWW.HAWSO.COM

ECN: 6210	REV. ECN: 6305	BY: JU	MODEL(S)	PART NUMBER
DRAWN: JU	DATE: 01/13/25	CHKD: JL	9006, PORTABLE EYEWASH STAND	0510002288.D
APPROVED: JU	DATE: 07/29/25	SCALE: 1:24	DRAWING TYPE: INSTALLATION	REVISION
			SIZE: A	2
				SHEET 2 OF 3

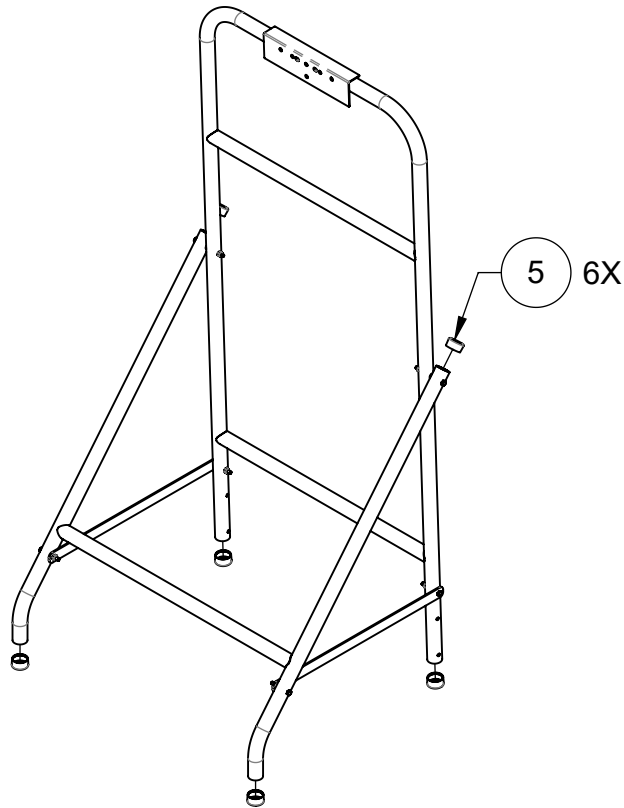

STEP 3

USE THE CAPS (ITEM 5) TO COVER THE OPEN TUBE ENDS.

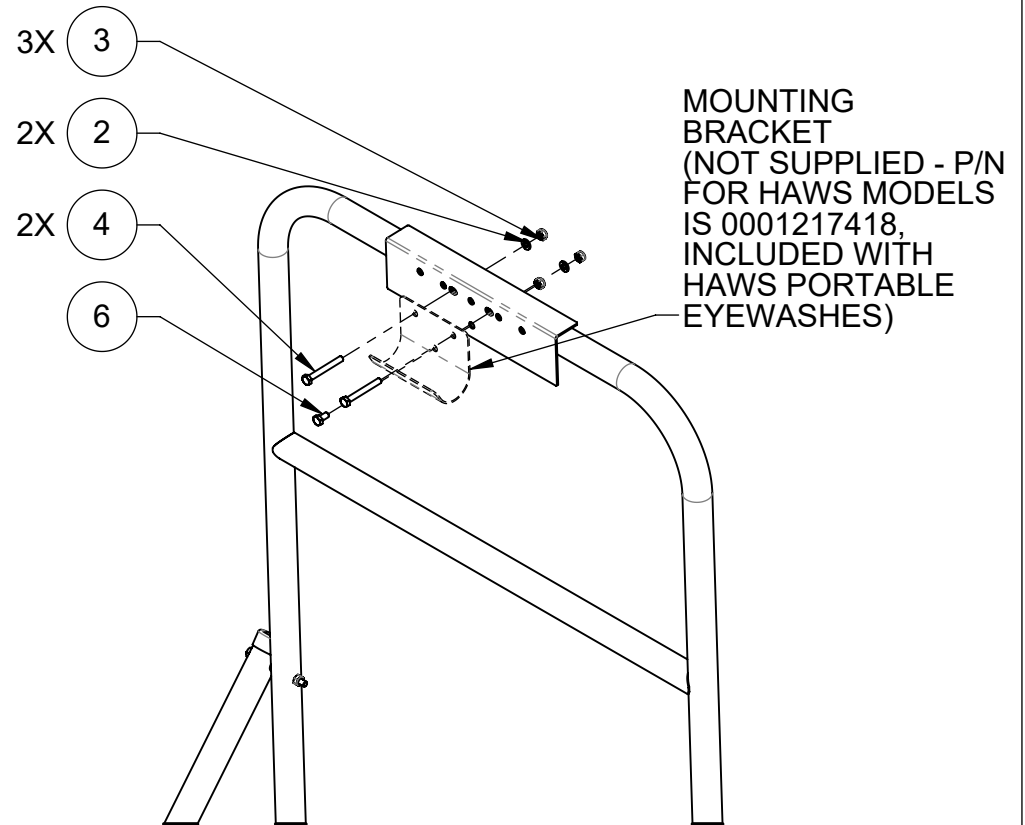

STEP 4

INSTALL THE PORTABLE EYEWASH MOUNTING BRACKET (NOT INCLUDED) USING MEDIUM SCREWS (ITEM 4), SHORT SCREW (ITEM 6), WASHERS (ITEM 2), AND NUTS (ITEM 3) AS NEEDED.



1455 KLEPPE LANE
SPARKS, NEVADA 89431
(775) 359-4712 FAX (775) 359-7424
E-MAIL: HAWS@HAWS.CO.COM
WEBSITE: WWW.HAWS.CO.COM

ECN: 6210	REV. ECN: 6305	BY: JU	MODEL(S) 9006, PORTABLE EYEWASH STAND	PART NUMBER 0510002288.D
DRAWN: JU	DATE: 01/13/25	CHKD: JL		REVISION 2
APPROVED: JU	DATE: 07/29/25	SCALE: 1:16	DRAWING TYPE: INSTALLATION	SIZE: A
				SHEET 3 OF 3